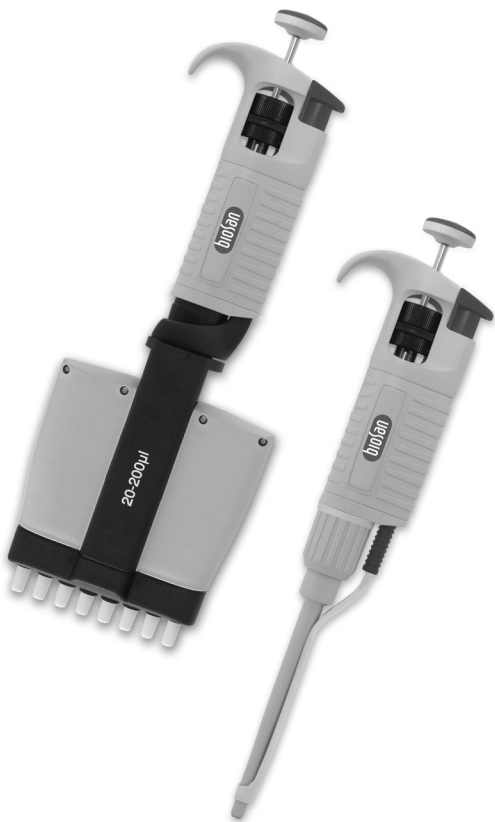
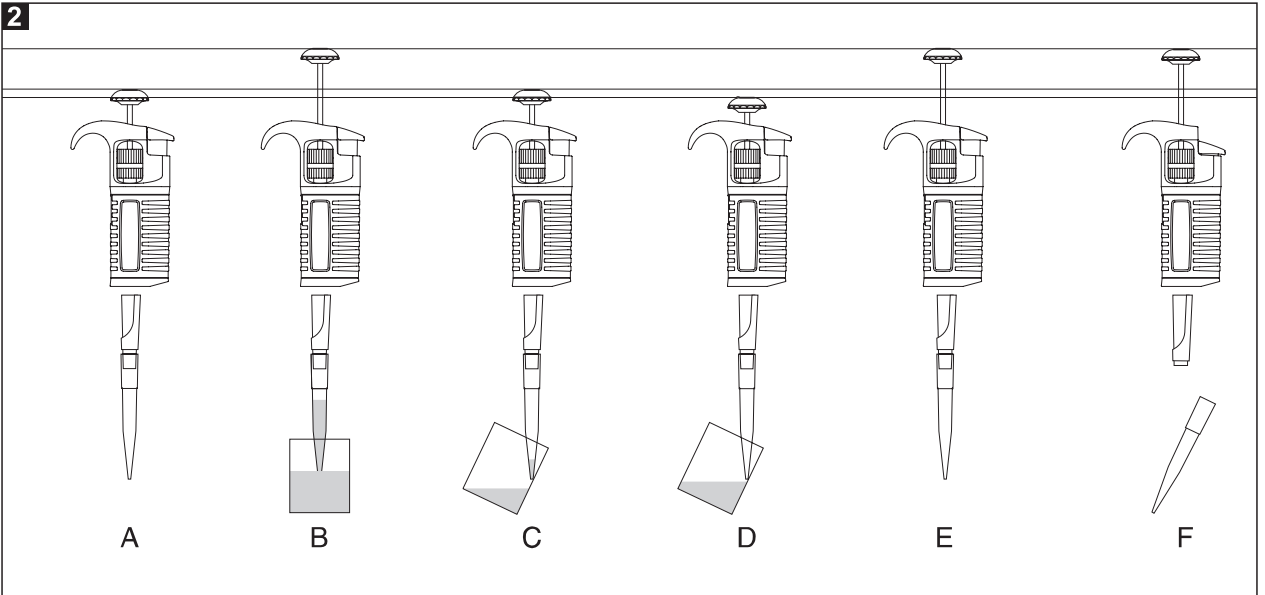
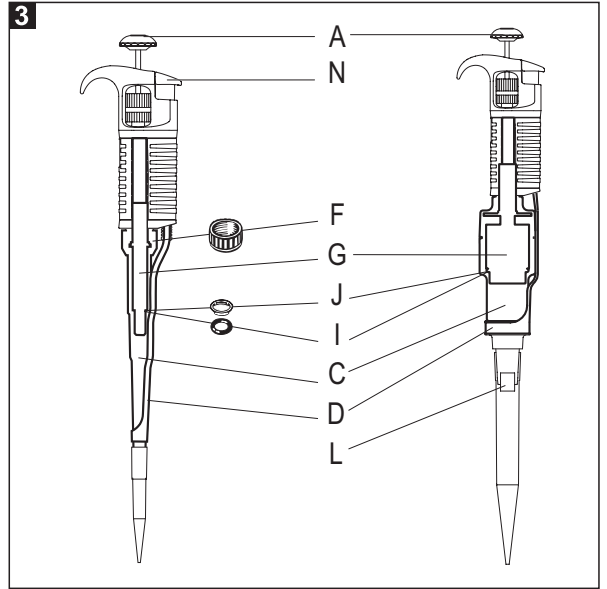
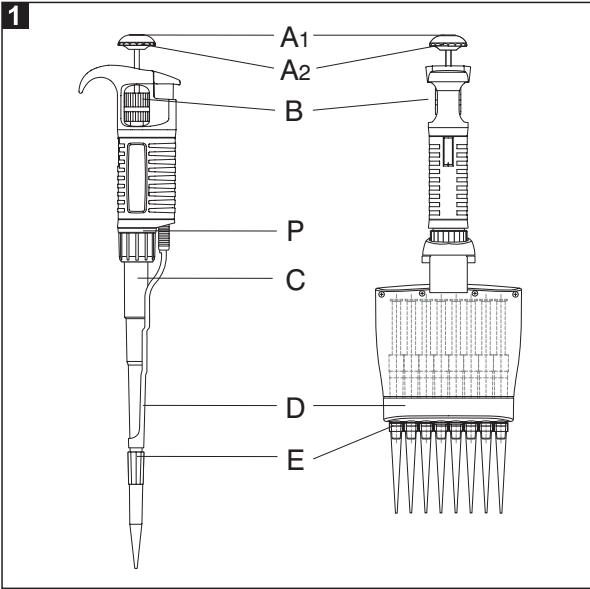
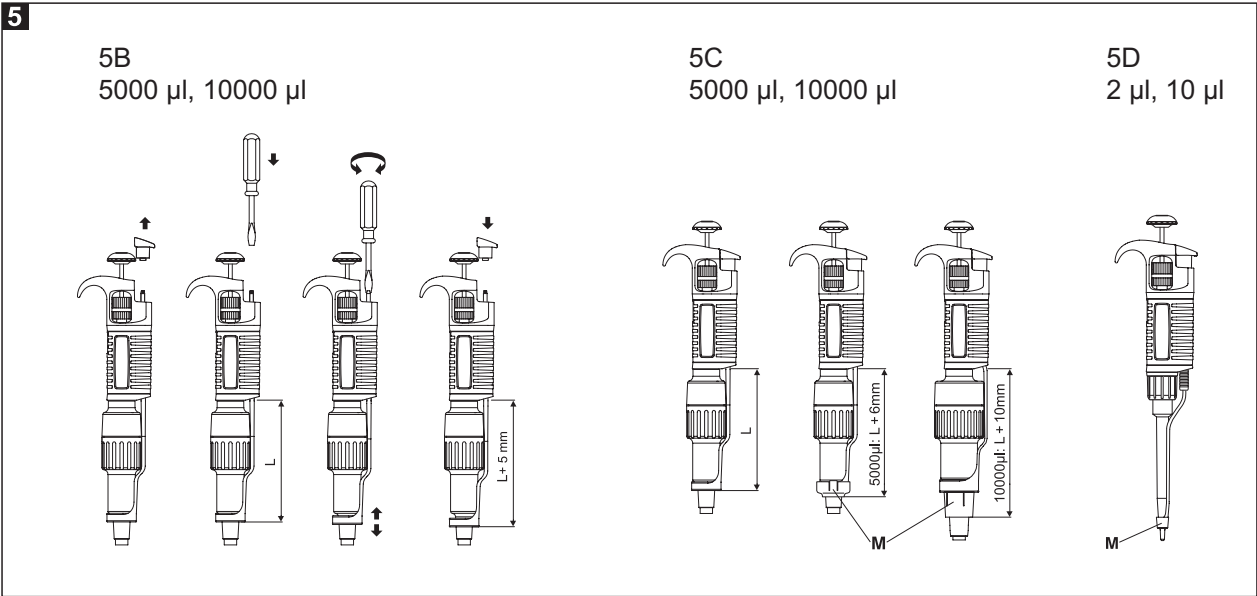
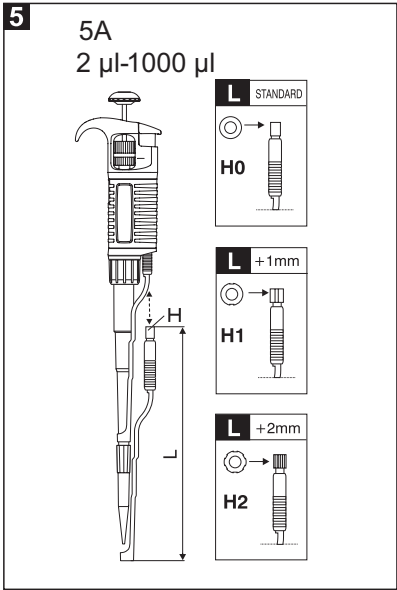
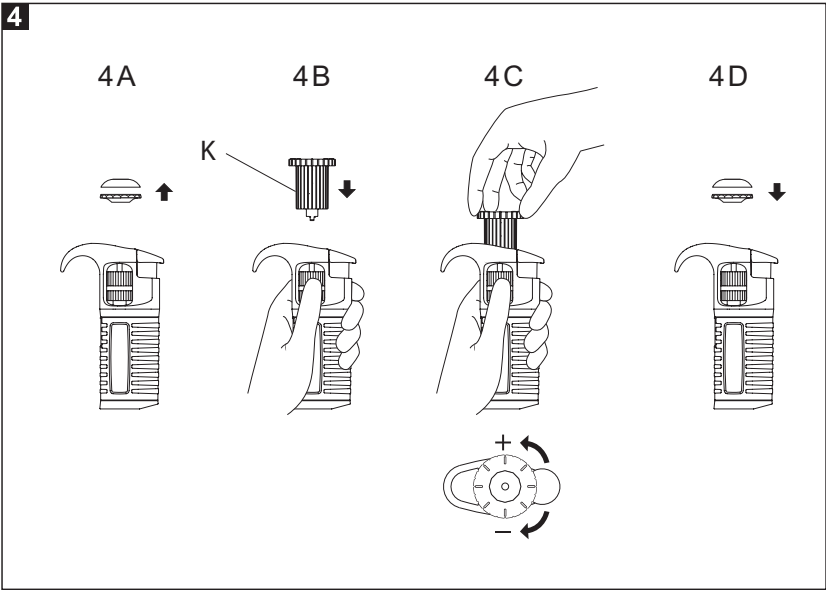
Assist series pipettes

ENGLISH	6 – 23
DEUTSCH	24 – 41
FRANÇAIS	42 – 59
ESPAÑOL	60 – 77
PORTUGUÊS	78 – 93
ITALIANO	94 – 109
РУССКИЙ	110 – 127

CONTENTS

- 1 – INTRODUCTION
- 2 – SETTING THE VOLUME
- 3 – ASPIRATING AND DISPENSING LIQUIDS
- 4 – PRE-RINSING
- 5 – DENSE AND VISCOUS LIQUIDS
- 6 – PIPETTE TIPS
- 7 – RECOMMENDATIONS
- 8 – RECALIBRATION
- 9 – TROUBLESHOOTING
- 10 – CLEANING AND STERILIZATION
- 11 – PIPETTE KIT AND ACCESSORIES
- 12 – SPARE PARTS

1 – INTRODUCTION

The **Assist** pipette is a volumetric instrument designed to measure and transfer liquids precisely and safely. It can accommodate volumes from 0.1 µl to 10,000 µl depending on the model.

The multichannel pipettes have been designed for the filling of laboratory microplates. The pipettes enable precise and simultaneous delivery of 8 or 12 preset-volume doses of liquid. The multichannel pipettes are produced in four ranges of volumes: 0,5-10 µl, 5-50 µl, 20-200 µl, 50-300 µl.

All **Assist** pipettes feature a analog counter which shows the pipetting volume. The set volume is visible in the window on the handle. Volumes are set by either turning the pipetting pushbutton (Fig. 1A2) or the adjustment knob (Fig. 1B) in the appropriate direction.

Liquids are aspirated into disposable tips attached to the pipette shaft. Disposable tips ensure maximum safety and eliminate the possibility of crosscontamination of the liquid samples.

To protect the user from contaminated tips, the **Assist** pipette is equipped with a built-in tip ejector, (Fig. 1D).

The construction of the ejector enables the user to set up the length. The adjustable tip ejector accommodates every variety of tips available on the market. When using narrow tubes, it may be necessary to remove the tip ejector. It is simply removed by pulling down.

Single channel pipette's ejector

Adjusting of the Tip Ejector Length

- in 2-1,000 µl pipettes (Fig. 5A):

The regulated "H" spacers, allow for regulating the length of tip ejector by +1 mm or +2 mm. An "H0" spacer is inserted on a standard basis.

- in 5,000 and 10,000 µl pipettes (Fig. 5B):

The length of tip ejector is adjusted by screwing in or screwing out the tip ejector rod with a small screwdriver. Turn the screwdriver counter clockwise to increase the length of tip ejector, and clockwise to reduce the length of tip ejector. The ejector can be adjusted by up to 5 mm.

If the ejector adjustment described above is not sufficient or the diameter of the ejector opening is too large to eject the tip it will be necessary to put the ejector cap "M" onto the ejector, (Fig. 5C).

- in 2 and 10 µl pipettes (Fig. 5D)

The 2 µl - 10 µl pipettes may require the use of an ejector cap „M” to efficiently eject certain brands of pipette tips. Simply place the cap, supplied with the pipette, on the bottom of the pipette shaft and slide the cap upwards until it surrounds the bottom of the tip ejector.

Multichannel pipette's ejector

The multichannel pipette construction allows to rotate the multichannel module and to eject the tips through 360°, therefore, the most convenient position may be selected to fill the microplates. It is recommended to turn the module clockwise. Due to the long stroke of the ejector in the multichannel pipette, the adjustment of its length is not necessary.

The **Assist** pipette is a high quality instrument which offers excellent accuracy and precision.

The accuracy and precision (repeatability) of liquid volume depends on the quality of tips used. The values for accuracy and precision shown in the table below were obtained using compatible tips. Only these tips guarantee correct operation of the pipettes and ensure accuracy and precision of liquid sampling.

The single channel pipettes

Model	Volume [µl]	Accuracy [%]	Precision [%]
AP2	0.2	± 12.0	± 6.0
	1.0	± 2.7	± 1.3
	Max 2.0	± 1.5	± 0.7
AP10	Min 0.5	± 4.0	± 2.8
	5.0	± 1.0	± 0.6
	Max 10.0	± 0.5	± 0.4
AP20	Min 2	± 3.0	± 1.5
	10	± 1.0	± 0.5
	Max 20	± 0.8	± 0.3
AP50	Min 5	± 2.5	± 2.0
	25	± 1.0	± 0.6
	Max 50	± 0.8	± 0.4
AP100	Min 10	± 1.6	± 0.80
	50	± 0.8	± 0.24
	Max 100	± 0.8	± 0.20
AP200	Min 20	± 1.2	± 0.60
	100	± 0.8	± 0.25
	Max 200	± 0.6	± 0.20
AP250	Min 50	± 1.0	± 0.4
	125	± 0.8	± 0.3
	Max 250	± 0.6	± 0.3

Model	Volume [µl]	Accuracy [%]	Precision [%]
AP1000	Min 100	± 1.6	± 0.40
	500	± 0.7	± 0.20
	Max 1,000	± 0.6	± 0.15
AP5000	Min 500	± 1.2	± 0.50
	2,500	± 0.6	± 0.20
	Max 5,000	± 0.5	± 0.15
AP10000	Min 1,000	± 2.5	± 0.6
	5000	± 0.8	± 0.3
	Max 10,000	± 0.5	± 0.2

The multichannel pipettes

Model	Volume [µl]	Accuracy [%]	Precision [%]
AP8-10 AP12-10	Min 0.5	±10.0	± 8.0
	Max 10	±4.0 ±2.0	± 2.0 ± 1.2
AP8-50 AP12-50	Min 5	±4.0	± 2.5
	Max 25 50	±3.0 ±1.6	± 1.2 ± 0.6
AP8-200 AP12-200	Min 20	±3.0	± 1.5
	Max 100 200	±1.5 ±1.0	± 0.8 ± 0.6
AP8-300 AP12-300	Min 50	±1.6	± 1.5
	Max 150 300	±1.2 ±1.0	± 1.0 ± 0.6

The accuracy and precision are obtained gravimetrically using compatible tips, performing at least 10 measurements of distilled water at a temperature of 20±1°C, according to EN ISO 8655 standard.

The design enables the user to recalibrate the pipette themselves according to the rules described in section 8.

2 – SETTING THE VOLUME

Setting the volume of the aspirated liquid can be done either using the pipetting pushbutton (Fig. 1A2) or by the adjustment knob (Fig. 1B). The volume display shown by the counter has three digits. These should be read from top to bottom. The lowest counter dial shows a scale which allows volume setting in the minimum graduation range. Examples of the black and red digits:

The single channel pipettes

Model	Counter Readings	Set Volume	Basic Degree			
AP2	<table border="1"><tr><td>1</td></tr><tr><td>2</td></tr><tr><td>5</td></tr></table>	1	2	5	1.25 µl	0.002 µl
1						
2						
5						
AP10	<table border="1"><tr><td>0</td></tr><tr><td>7</td></tr><tr><td>5</td></tr></table>	0	7	5	7.5 µl	0.02 µl
0						
7						
5						
AP20	<table border="1"><tr><td>1</td></tr><tr><td>2</td></tr><tr><td>5</td></tr></table>	1	2	5	12.5 µl	0.02 µl
1						
2						
5						
AP100	<table border="1"><tr><td>0</td></tr><tr><td>7</td></tr><tr><td>5</td></tr></table>	0	7	5	75 µl	0.2 µl
0						
7						
5						
AP200	<table border="1"><tr><td>1</td></tr><tr><td>2</td></tr><tr><td>5</td></tr></table>	1	2	5	125 µl	0.2 µl
1						
2						
5						
AP1000	<table border="1"><tr><td>0</td></tr><tr><td>7</td></tr><tr><td>5</td></tr></table>	0	7	5	0.75 ml	2 µl
0						
7						
5						
AP5000	<table border="1"><tr><td>1</td></tr><tr><td>2</td></tr><tr><td>5</td></tr></table>	1	2	5	1.25 ml	10 µl
1						
2						
5						
AP10000	<table border="1"><tr><td>0</td></tr><tr><td>7</td></tr><tr><td>5</td></tr></table>	0	7	5	7.5 ml	20 µl
0						
7						
5						

The multichannel pipettes

Model	Counter Readings	Set Volume	Basic Degree			
AP8-10 AP12-10	<table border="1"><tr><td>0</td></tr><tr><td>3</td></tr><tr><td>5</td></tr></table>	0	3	5	3.5 µl	0.02 µl
0						
3						
5						
AP8-50 AP12-50	<table border="1"><tr><td>0</td></tr><tr><td>6</td></tr><tr><td>5</td></tr></table>	0	6	5	6.5 µl	0.10 µl
0						
6						
5						
AP8-200 AP12-200	<table border="1"><tr><td>0</td></tr><tr><td>8</td></tr><tr><td>5</td></tr></table>	0	8	5	85.0 µl	0.20 µl
0						
8						
5						
AP8-300 AP12-300	<table border="1"><tr><td>2</td></tr><tr><td>5</td></tr><tr><td>0</td></tr></table>	2	5	0	250.0 µl	1.0 µl
2						
5						
0						

To attain the maximum accuracy, set volume must be approached from a higher value.

If the desired value is lower than the previous value, simply adjust the value down to 1/3 turn above the required setting.

If the desired value is higher than the previous value, adjust the value to 1/3 turn above the required value. The setting should then be performed in the following manner:

Slowly turn the pipetting pushbutton knob or the black adjustment knob to the required volume.

If the knob is accidentally turned too far, below the required value, the process must be repeated.

3 – ASPIRATING AND DISPENSING LIQUIDS

The single channel pipettes

Place a tip on the shaft of the pipette. See Section 6 for the appropriate tip. Press the tip on firmly using a slight twisting motion. This will ensure an airtight seal.

The multichannel pipettes

In order to put tips on the multichannel shafts hold the pipette vertically and press it against the tips in the rack box, until the shafts retract about 1.5 mm into the multichannel module. The suspension system of the shafts ensures even and tight sealing of tips. The rocking movement does not have to be performed to seal the tips tightly.

Important: Never aspirate liquids into the Assist pipette without a tip attached.

Aspiration

Press the pushbutton to the first positive stop, (Fig. 2A). Holding the pipette vertically, immerse the tip into the sample liquid. The depth to which the tip is immersed in the sample liquid depends on the model.

Model	Immersion depth [mm]
AP2	≤ 1
AP10	≤ 1
AP20, AP50, AP100	2 - 3
AP200, AP250, AP1000	2 - 3
AP5000	3 - 6
AP10000	5 - 7

Release the pushbutton slowly and smoothly to aspirate the sample, (Fig. 2B). Wait one second and then withdraw the tip from the liquid. If the pipette tip is not immersed to the recommended depth or if the pipetting pushbutton is rapidly released, air may enter the disposable tip.

Avoid touching the opening of the tip.

Dispensing:

- Place the end of the tip against the inside wall of the vessel at an angle of 10° to 40°.
- Press the pushbutton smoothly to the first stop, (Fig. 2C). Wait one second.
- Press the pushbutton to the second stop to expel any remaining liquid, (Fig. 2D).
- Keeping the pushbutton depressed to the very end, remove the pipette by drawing the tip against the inside surface of the receiving vessel.
- Release the pushbutton to its starting position, (Fig. 2E).
- Eject the tip by pressing the tip ejector button, (Fig. 2F). Remember to change the tip whenever a different kind of liquid is to be sampled.

Filters

A replaceable filter, fitted in a seat in the bottom part of the shaft, is used in 5,000 µl and 10,000 µl pipettes (Fig. 3L). This filter prevents the aspirated liquid from entering into the shaft and thus from contaminating the inside of the shaft and the piston. Using the filter is particularly important when aspirating and dispensing large volumes of liquid.

If the filter becomes wet during liquid aspiration it should be replaced with a new one.

4 – PRE-RINSING

When pipetting liquids of higher viscosity or lower surface tension than water (e.g. sera or organic solvents), a film of liquid is formed on the inside wall of the pipette tip. This film can create an error. As the film remains relatively constant in successive pipetting operations with the same tip, this error can be eliminated by forming the film before transferring the first sample. This is done by aspirating a sample and dispensing it back into the same vessel. Now, as the film is already formed, all of the following samples will have better accuracy and repeatability.

This pre-rinsing operation should be repeated when the volume to be aspirated is changed or when a new tip is used.

5 – DENSE AND VISCOUS LIQUIDS

Assist pipette specifications of accuracy and precision are based on pipetting distilled water. Handling of liquids with physical qualities of density, viscosity and surface tension

which differ substantially from water may need gravimetrically checked compensation of the volume setting. Normally the degree of error resulting from heavy or viscous liquids is negligible if pipetting is done slowly and carefully. It is most important to give the liquids some time to react to the change of pressure by holding the pipette tip in its position for at least 2 sec. after the aspiration and the blow out stroke.

If in extreme cases, this method of operation does not result in accurate values, a compensation may be achieved as follows: Weigh the liquid pipetted when the **Assist** pipette is set to the nominal value. Then calculate the set-off from the nominal value:

$$\text{Correction, value} = 2 \times \text{nominal val.} \frac{m}{\gamma}$$

m – weight of the sample
 γ – density of liquid

Check this operation once again and correct if necessary. Note the corrected value for further pipetting of the same kind of liquid.

6 – PIPETTE TIPS

Our recommended tips are made from high performance polypropylene and their quality guarantees the precision and accuracy associated with the **Assist** pipette. Strict control is maintained throughout the manufacturing process to ensure the highest quality.

The accuracy and precision figures for the **Assist** pipette are only guaranteed when our recommended tips are used. The use of inferior quality tips will seriously degrade the performance of the **Assist** pipette.

Tips 10

These tips are used for volumes between 0.1 µl and 10 µl. They are used with the AP2, AP10, AP8-10 & AP12-10 models.

Tips 200

These tips are used for volumes between 2 µl and 200 µl. They are used with the AP20, AP50, AP100, AP200, AP8-50, AP12-50, AP8-200 and AP12-200 models.

Tips 300

These tips are used for volumes between 50 µl and 300 µl. They are used with the AP250, AP8-300 and AP12-300 models.

Tips 1,000

These tips are used for volumes between 100 µl and 1,000 µl. They are used with the AP1000 model.

Tips 5,000

These tips are used for volumes between 500 µl and 5,000 µl. They are used with the AP5000 model.

Tips 10000

These tips are used for volumes between 1,000 and 10,000 µl. They are used with the AP10000 model.

7 - RECOMMENDATIONS

Observing the following recommendations will ensure maximum possible accuracy and precision of liquid sampling.

- Make sure to operate the **Assist** pipette slowly and smoothly.
- The depth of immersion in the sample liquid should be the minimum necessary and should remain constant during aspiration.
- The **Assist** pipette should be held in a vertical position.
- Change the tip when volume setting is changed or when a different liquid is to be aspirated.
- Change the tip if a droplet remains on the end of the tip from the previous pipetting operation.
- Each new tip should be pre-rinsed with the liquid to be pipetted.
- Liquid should never enter the **Assist** pipette shaft. To prevent this:
 - Press and release the pushbutton slowly and smoothly.
 - Never turn the pipette upside down.
 - Never lay the pipette on its side when there is liquid in the tip.
- Never force the volume setting beyond its recommended limits.
- When pipetting liquids with temperatures different from the ambient temperature, it is recommended to pre-rinse the tip several times before use.
- Do not pipette liquids with temperatures above 70°C.
- When pipetting acids or corrosive solutions which emit vapours, it is recommended to disassemble the shaft and rinse the piston and seal with distilled water after finishing the pipetting operation.

8 – RECALIBRATION

Pipettes are calibrated by gravimetrically, using our recommended tips and distilled water, at a temperature of 20±1°C, according to EN ISO 8655 standard.

If during pipette operation you find that the accuracy error (the difference between the real aspirated volume and the preset volume) exceeds the permissible value given in the table in section 1, the pipette recalibration procedure should be carried out.

Before starting the recalibration it is necessary to check whether the following requirements have been met during error determination:

- the ambient temperature, and the temperature of the pipette, tips and water was identical
- the density of the liquid used is close to that of distilled water
- a balance with appropriate sensitivity has been used.

Volume checked [µl]	Balance sensitivity [mg]
0.1 - 10	≤0.001
10 - 100	≤0.01
>100	≤0.1

- mg/µl conversion factor has been taken into account
 - the requirements shown in sections 3 and 7 have been met
- If the above conditions are satisfied and the accuracy error for selected volume given in section 1 exceeds the permissible value, the pipette recalibration procedure should be carried out.

The recalibration can be performed within one full turn of the key to the right or to the left only.

Recalibration conditions:

Ambient temperature, the temperature of the pipette, tips and liquid should be within the range 20-25°C and stabilised during weighing within ±0.5°C

- Measurements should be conducted using distilled water
- Balance sensitivity should be suitable for the volume to be measured

Recalibration procedure:

- Set the dose volume depending on the pipette volume according to the following table:

The single channel pipettes

Model	Range of the pipette volumes [μl]	Preset volume [μl]	Permissible volumes [μl]	Volume change
				ΔV for full turn of the calibration key [μl] (24 increments)
AP2	0.1 - 2	0.2	0.176 - 0.224	0.06
AP10	0.5 - 10	0.5	0.48 - 0.52	0.33
AP20	2 - 20	2	1.94 - 2.06	0.63
AP50	5 - 50	5	4.875 - 5.125	2.50
AP100	10 - 100	10	9.84 - 10.16	2.50
AP200	20 - 200	20	19.76 - 20.24	6.30
AP250	50 - 250	50	49.5 - 50.5	6.30
AP1000	100 - 1000	100	98.4 - 101.6	25.00
AP5000	500 - 5,000	500	494 - 506	125.00
AP10000	1,000 - 10,000	1,000	975 - 1,025	250.00

The multichannel pipettes

Model	Range of the pipette volumes [μl]	Preset volume [μl]	Permissible volumes [μl]	Volume change
				ΔV for full turn of the calibration key [μl] (24 increments)
AP8-10 AP12-10	0.5 - 10	0.5	0.45 - 0.55	0.33
AP8-50 AP12-50	5 - 50	5	4.8 - 5.2	1.67
AP8-200 AP12-200	20 - 200	20	19.4 - 20.6	6.30
AP8-300 AP12-300	50 - 300	50	49.2 - 50.8	10.00

- Perform 5 aspirations, weigh each one and calculate the average value of the aspirations
- Perform three aspiration series (each series should include the aspirations from all channels), weigh each time and calculate the average value of the aspirations.
- Calculate average aspirated volume in μl multiplying the average aspiration amount [mg] by the distilled water density coefficient [μl/mg], which depends on temperature and pressure according to the following table:

Temperature [°C]	Pressure [kPa]		
	95.0	101.3	105.0
20	1.0028	1.0029	1.0029
21	1.0030	1.0031	1.0031
22	1.0032	1.0033	1.0033
23	1.0034	1.0035	1.0036
24	1.0037	1.0038	1.0038
25	1.0039	1.0040	1.0040

If the average aspirated volume exceeds the permissible value, the following should be done:

- Remove the pipetting pushbutton, (Fig. 4A),

Warning: The pipetting pushbutton consists of 2 parts: the lower volume-setting section (Fig. 1A2) and the upperbutton (Fig. 1A1). After removal of the pushbutton, both parts will separate.

- Holding the volume setting knob to protect it against rotation, insert the calibration key into the cuts of the calibration screw, (Fig. 4B),
- Turn the key clockwise to reduce the aspirated volume, or counter-clockwise to increase the volume. One full turn of the calibration key changes the pipette aspiration volume by the amount given in the table, (Fig. 4C),
- Take out the key and fix the pipetting pushbutton (Fig. 4D). The pipetting pushbutton should be fixed by placing the lower section on the arbor first (Fig. 1A2) and then the button (Fig. 1A1).

Determine the average aspirated volume. The average volume should be within the permissible range shown in the table. If the volume exceeds the values stated, the recalibration procedure should be repeated.

When pipetting liquids with physical properties considerably different from those of water, follow the rules as described in section 5.

9 – TROUBLESHOOTING

If you notice an improper pipette operation identify the cause and eliminate the fault. To do this, follow the instruction accordingly. Replacement of parts should be required only occasionally, and should not occur under normal pipette use.

Problem	Cause	Solution
Droplets of liquid remain in the pipette tip.	The pipette tip is emptied too fast.	Decrease the speed of pressing the pipette pushbutton.
	The tip wettability has increased due to extensive use.	Replace the tip with a new one.
Droplets of air appear in the liquid aspirated into the tip.	The pipette tip immersion is too shallow.	Immerse the tip deeper according to the instructions.
	The pipette tip is incorrectly pressed onto the pipette shaft.	Press the pipette tip firmly.
	The tip is damaged or worn out due to extensive use.	Replace the tip with a new one.
The pipette incorrectly aspirates the liquid or liquid drops out from the tip.	The pipette tip is incorrectly pressed onto the pipette shaft.	Press the pipette tip firmly.
	The shaft nut is loose (Fig. 3F) in the models 2 - 1000	Tighten the shaft nut.
	The sealing surface of the shaft is cracked or scored.	Remove the tip ejector. Unscrew the shaft nut, inspect the shaft and the piston assembly. Replace the damaged parts (see Section 12). When reassembling the pipette, the nut should be hand tightened. In the models AP2, AP10 and AP20, the damage of the shaft may also cause a damage of the piston assembly. Replace the damaged parts (see Section 12). When reassembling the pipette, the nut should be hand tightened. To remove the tip ejector in models AP5000 and AP10000, remove the ejector pushbutton (Fig. 3N) and using a screwdriver unscrew the tip ejector by turning the screwdriver counter-clockwise.

Problem	Cause	Solution
The pipette incorrectly aspirates the liquid or liquid drops out from the tip.	Damage to the piston or seal due to prolonged use with the aggressive liquids.	Disassemble the pipette as described above. Replace the piston, seal and O-ring (see Section 12). Rinse the inside of the shaft in distilled water and dry. Lubricate the seal and O-ring with the lubricant, that is included with each pipette. Following piston replacement, a calibration procedure is required. Note: The parts of AP2 and AP10 pipette should be lubricated evenly with a minimum amount of lubricant.
	The pipette is reassembled improperly.	Disassemble the pipette and reassemble it, observing the proper sequence of steps (Fig. 3).
	No lubricant on the sealing elements.	Remove the tip ejector. Unscrew the shaft nut, remove the shaft, piston assembly, seal and O-ring. Rinse the removed parts in distilled water and dry thoroughly. Lightly lubricate the inside surfaces of the seal and the O-ring with the included lubricant. Reassemble the pipette in the reverse order.
	Contamination of the inside of the pipette caused by extensive aspiration of chemically aggressive liquids or because liquid got inside the pipette.	Remove the tip ejector. Unscrew the nut, remove the shaft, piston assembly, seal and O-ring. Rinse the removed parts with distilled water and dry thoroughly. Lightly lubricate the inside surfaces of the seal and the O-ring with the lubricant. Reassemble the pipette in the reverse order.

If you find an increase in the pipetting force, which could happen after repetitive autoclaving of the pipette:

Remove the ejector. Unscrew the shaft nut, and then remove the shaft, piston assembly, seal and O-ring. Rinse the removed parts in distilled water and dry. Lubricate the internal surfaces of the seal and O-ring with lubricant that has been included with each pipette. Reassemble the pipette in opposite order.

Note: All parts of the pipette can be autoclaved (see Section 10). The shaft of the AP5000 and AP10000 models should be autoclaved without the filter.

If the problem continues after carrying out the above steps, contact your **Assist** pipette representative.

Before returning the pipette, please ensure that the pipette is completely free of any chemical, radioactive or microbiological contamination which could pose a threat during transport and repair. As far as it is possible, clean the pipette.

10 – CLEANING AND STERILISATION

Cleaning:

External surfaces of the pipetting pushbutton, the ejector pushbutton, the handgrip, the shaft nut and the adjustment knob may be cleaned using a cloth dampened in isopropyl alcohol. The remaining parts removed from the pipette during pipette disassembly may be washed with distilled water or isopropyl alcohol.

Warning:

Before using cleaning agents other than those recommended by the manufacturer, check the compatibility charts and consider chemical resistance of the following plastics: PP, PC, POM, PA, PPS, PVDF used to produce the pipette parts.

Sterilization:

The pipette can be sterilised in an autoclave at 121°C for 20 minutes. After sterilisation, the pipette should be dried and cooled to room temperature

It is recommended:

- to sterilize the pipettes in autoclave with an initial vacuum and drying cycle,
- prior to sterilization unscrew the shaft nut slightly in the AP2 to AP1000 pipettes, and unscrew the shaft slightly in the AP5000 and AP10000. After autoclaving these parts should be screwed tight again.
- prior to sterilization the multichannel pipette unscrew the nut connecting the handle and module slightly. After autoclaving these parts should be screwed tight again.

The shaft of the 5000 and 10000 models should be autoclaved without the filter.

The precision of the results should not alter if the pipetting process and autoclaving are carried out as described in this manual.

Because a slight change in the accuracy of the dosage may occur, it is recommended to:

- check the calibration of the pipette after the initial first, third and fifth autoclaving cycles and then after every 10 autoclaving cycles.

The pipettes are UV resistant which was confirmed by our tests. The recommended distance from the radiation source to exposed element should be not less than 50 cm. Prolonged and very intense UV exposure can cause de-coloration of pipette parts, without affecting its performance.

11 – PIPETTE KIT AND ACCESSORIES

Pipette kit:

The pipettes are delivered in the kits including:

- Pipette
- Instruction manual
- Calibration key
- Ejector length adjustment spacers (for pipette models AP2 to AP1000)
- Ejector cap (for pipette models AP2, AP10, AP5000, AP10000)
- Filters (for pipette models AP5000, AP10000)
- Lubricant
- Color Identification Rings

Accessories:

Description	Cat. no.
Carousel stand	BS-010522
Multiple stand	BS-010523
1-position stand	BS-010524
4-position stand	BS-010525

12 – SPARE PARTS

The spare parts indicated in Fig. 3, 4 and 5 that is:

- A: Pipetting pushbutton A1: Pushbutton A2: Knob
- B: Adjustment knob
- C: Shaft
- D: Ejector
- F: Shaft nut
- G: Piston assembly
- H: Spacer

- I: O-ring
- J: Seal
- K: Calibration key
- L: Filter
- M: Ejector cap
- N: Ejector pushbutton
- P: Identification ring

can be ordered from your **BIOSAN** representative (type of pipette and name of the part for this pipette should be specified).

To find your nearest service center and for technical support, visit <http://www.biosan.lv>

Warning: The replacement of the piston requires a calibration procedure to be carried out, as described in section 8.

All rights reserved. Product described in this manual is subject to availability and technical modification. Errors excepted.
 BIOSAN reserves the right to improve, enhance or otherwise modify its products without prior notification.

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